

# Instructions for downloading the NMR spectra from a network server for Windows 10 users - creating a new network location

1. In the lower left corner of your screen, in the search area type „map" and select the offered option „map network drive"

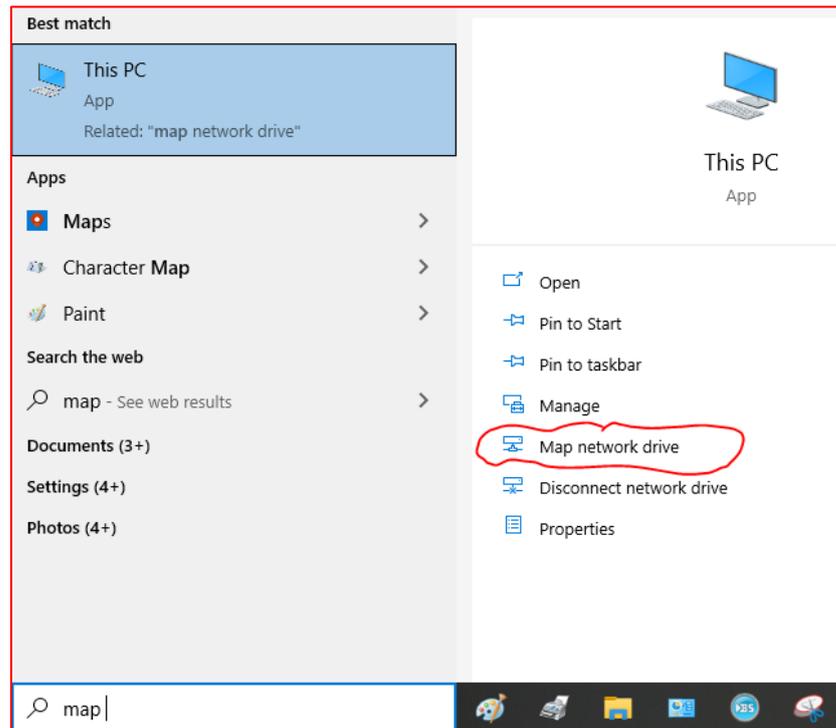


Figure 1.

2. When the window opens, select the option „Connect to a web site that you can use to store your documents and pictures“

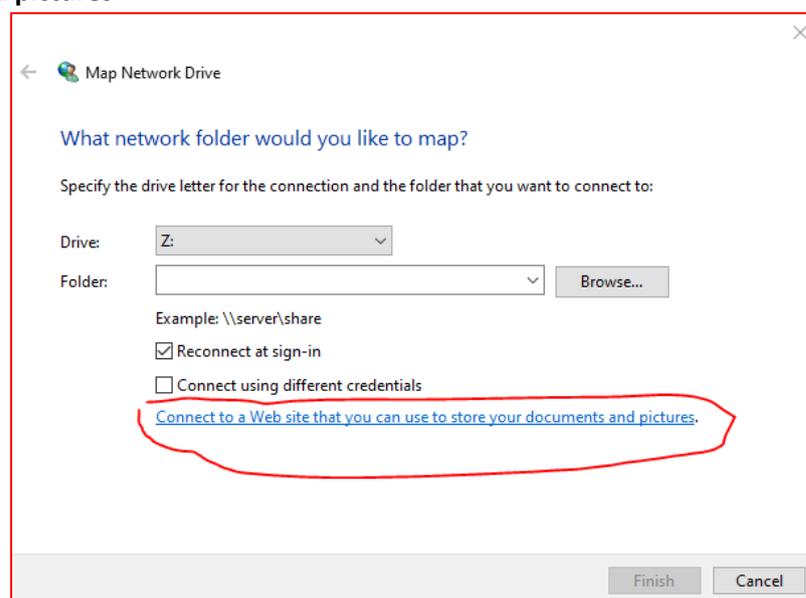


Figure 2.

3. Press „**Next**“ **three times** until you reach the window where you will enter the server address in the form: <https://nmr.irb.hr/username>

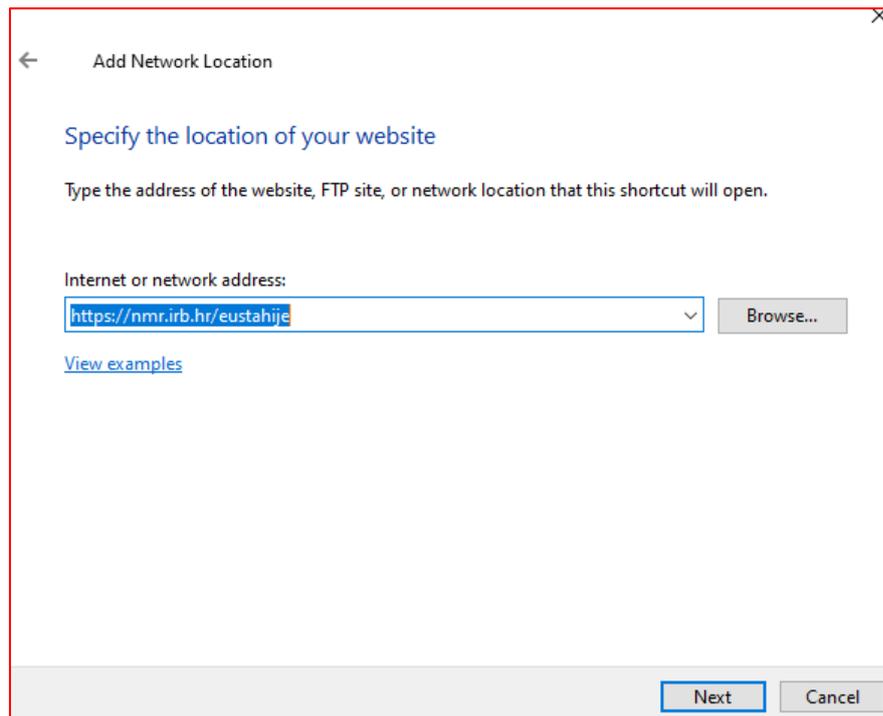


Figure 3.

4. After clicking „**Next**“, the system will ask for the password:


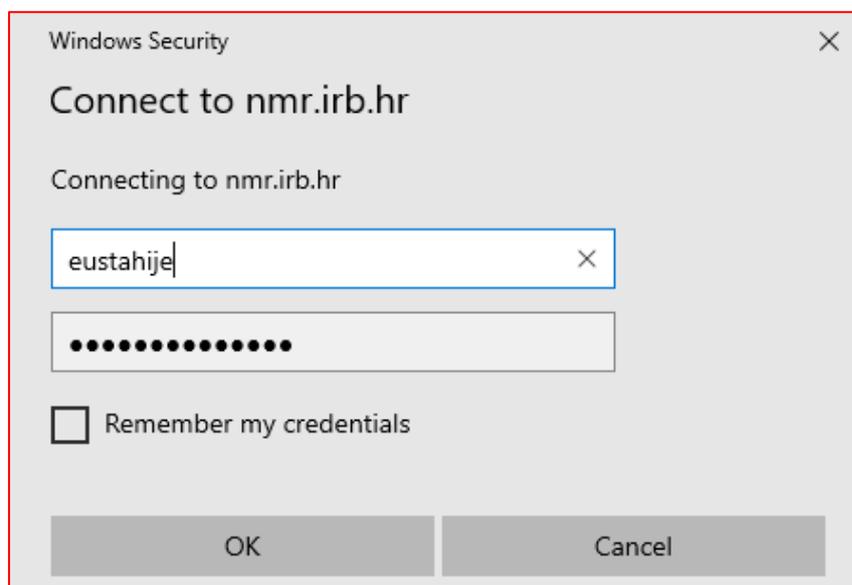


Figure 4.

5. Enter your password and press „**OK**“


- In the next step, the system will ask you to create a name for the new location. After selecting and entering the name, click „**Next**“:

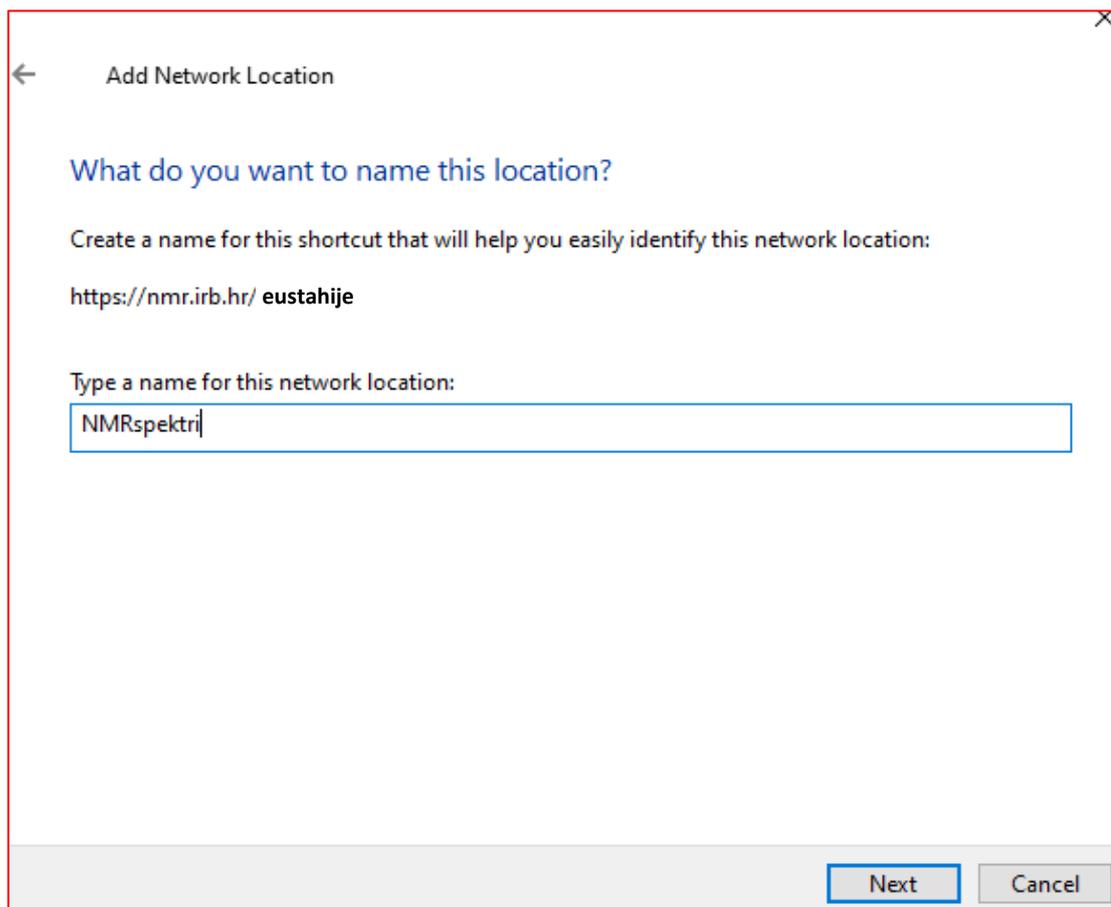

The screenshot shows a window titled "Add Network Location" with a back arrow on the left and a close button (X) on the right. The main text asks "What do you want to name this location?" and provides instructions: "Create a name for this shortcut that will help you easily identify this network location:" followed by the URL "https://nmr.irb.hr/ eustahije". Below this is a text input field with the label "Type a name for this network location:" and the text "NMRspektri" entered. At the bottom right, there are two buttons: "Next" (highlighted with a blue border) and "Cancel".

Figure 5.

- If you have successfully created a new network location, you will receive a notification:

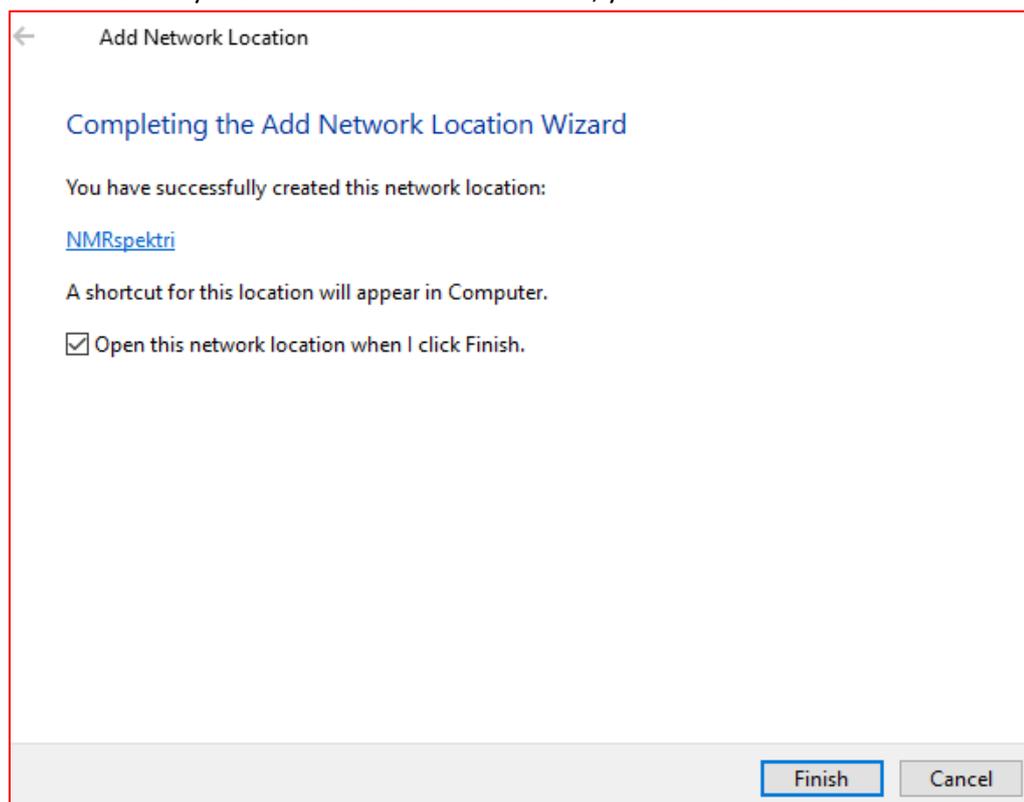

The screenshot shows a window titled "Add Network Location" with a back arrow on the left and a close button (X) on the right. The main text says "Completing the Add Network Location Wizard" and "You have successfully created this network location:" followed by the name "NMRspektri" in blue. Below this is the text "A shortcut for this location will appear in Computer." and a checked checkbox with the label "Open this network location when I click Finish." At the bottom right, there are two buttons: "Finish" (highlighted with a blue border) and "Cancel".

Figure 6.

8. Click „**Finish**” and the newly created network location will open making NMR spectra visible.
9. The new location will be visible each time you open Windows Explorer and select „**This PC**” as shown in Figure 7:

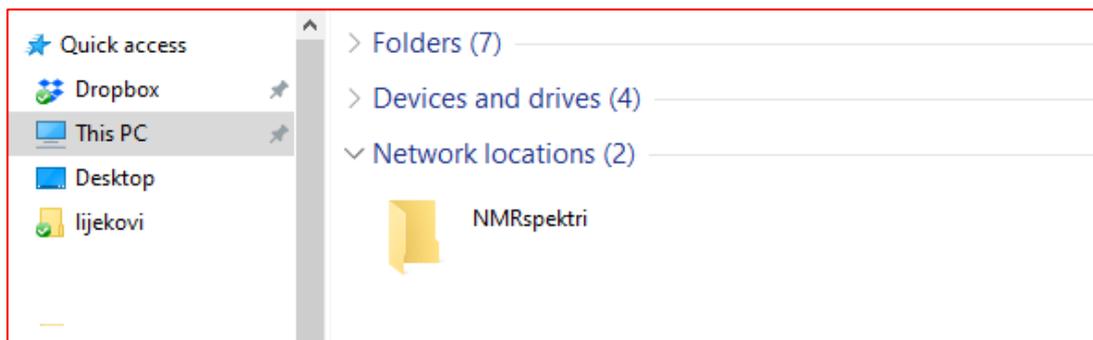


Figure 7.

From this location, you can copy spectra just like from any other location on your computer.

Please use your own username instead of fictional „eustahije” used here as an example. 😊